opensignals

(r)evolution

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CHANNEL TYPES

EMG	Electromyography
EDA	Electrodermal Activity
ECG	Electrocardiography
ACC	Accelerometer
LUX	Photodiode
RAW	Non-scaled signal visualization

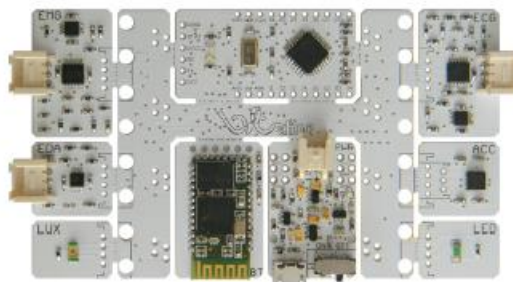
Fig. 17 BITalino Board

CHANNEL CONFIGURATION

BITalino's standard channel configuration:

- 1 EMG
- 2 EDA
- 3 ECG
- 4 ACC
- 5 LUX
- 6 Free for other (third party) components

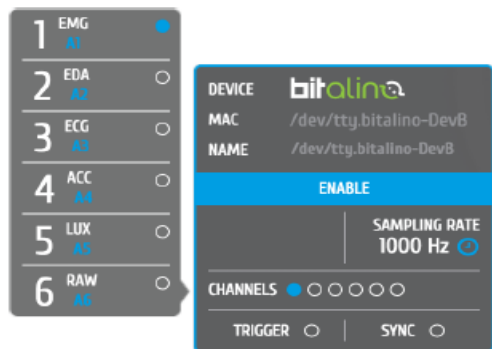
Fig. 18 BITalino extended Open Signals header

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LED STATES

White LED	Low frequency blinking on stand-by mode
White LED	High frequency blinking when acquiring
Red LED (MCU)	Battery is running low
Red LED (Power Unit)	BITalino is turned off; Battery's charging

5.3 BioSignalsPlux

CHANNEL TYPES:

ECC	Electrocardiography*
EMG	Electromyography*
HANDGR	Handgrip (Force Measurement)*
RESP	Respiration*
RAW	Non-scaled signal visualization
XYZ	Accelerometer
CUSTOM	User-scaled visualization
GONIO	Goniometer*
EDA	Electrodermal Activity*
BVP	Blood Volume Pressure*
EEG	Electroencephalography*

Fig. 19 8 channel BiosignalsPlux hub

*These channel types have pre-set signal visualization scales

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LED STATES (for single-device and multi-device acquisition)

Main Rules:

1 blink	Device is idle
2 blinks	Device is acquiring
Green Blink	Bluetooth on
Red Blink	Low Battery
Green	Charging
Orange blink	Internal acquisition

Detailed Information:

none	Power down (RTC is running)
1 Green Blink	Idle; Bluetooth on
2 Green Blinks	Real-time acquisition; Bluetooth on
1 Red Blink	Idle; Low battery; Bluetooth on
2 Red Blinks	Real-time acquisition; Low battery; Bluetooth on
1 Black Blink in the Background	Idle; Charging; Bluetooth on
2 Black Blinks in the Background	Real-time acquisition; Charging; Bluetooth on
1 Orange Blink	Waiting for external trigger to start internal acquisition; Bluetooth off

1 Orange Blink alternating with 1 Red Blink	Waiting for external trigger to start internal acquisition; Low battery; Bluetooth off
2 Orange Blinks	Internal acquisition; Bluetooth off
Orange - Red Blinks	Internal acquisition; Charging; Bluetooth off
2 Orange blinks in green Background	Internal acquisition; Charging; Bluetooth off
Orange - Green Blinks	Internal acquisition; Bluetooth on
Green - Red blinks	Internal acquisition; Low Battery; Bluetooth on
Orange - Black blinks in green background	Internal acquisition; Charging; Bluetooth on

BUTTON BEHAVIOUR

Device state before button press:	Device state after button press:
Power Down	Idle
Idle or real-time acquisition	Power down
Internal acquisition (Bluetooth on)	Intern acquisition (Bluetooth off)
Internal acquisition (Bluetooth off)	Internal Acquisition (Bluetooth on)

5.2.1 BiosignalsPlux Explorer Channels - 4 Channel Device

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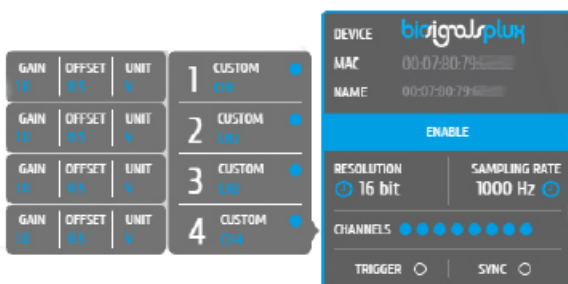
Fig. 20 4 channel BiosignalsPlux extended Open Signals header

5.2.2 BiosignalsPlux Researcher/Professional Channels - 8 Channel Device

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5.3 Motion Plux

CHANNEL TYPES:

ACC Accelerometer
MAG Magnetometer
RAW Non-scaled signal visualization
CUSTOM User scaled signal visualization

CHANNEL CONFIGURATION

MotionPlux' standard configuration:

- 1 Accelerometer (X-Axis)
- 2 Accelerometer (Y-Axis)
- 3 Accelerometer (Z-Axis)
- 4 Magnetometer (X-Axis)
- 5 Magnetometer (Y-Axis)
- 6 Magnetometer (Z-Axis)

Fig. 22 MotionPlux

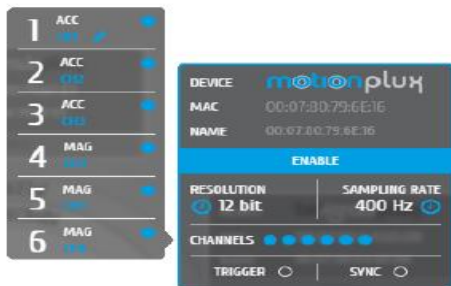
Fig. 23 MotionPlux extended Open Signals header

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LED STATES

Blinks 1/sec	Stand-By
Blinks 2/sec	Acquiring
Blinks Green	High battery level
Blinks Red	Low battery level
Constantly Green	Charging

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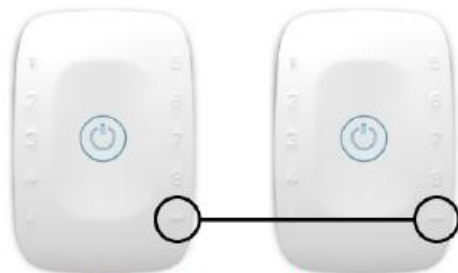
Fig. 24 Channels for sync cable connection

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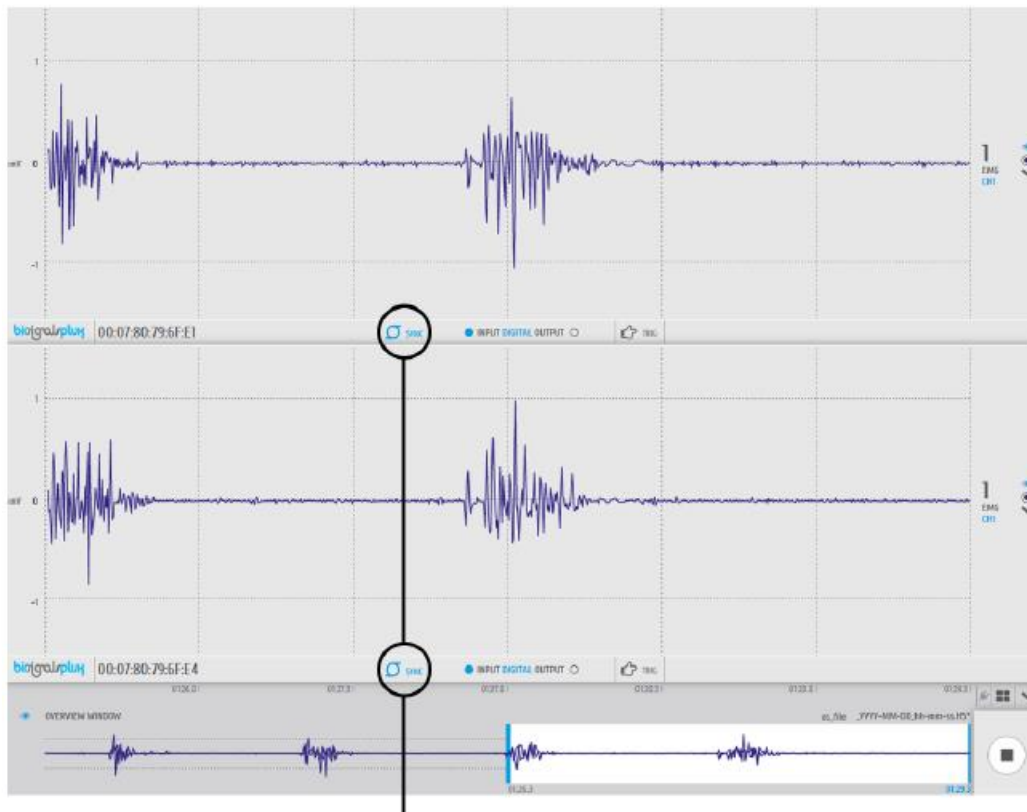
Example: Device A 5 samples in $\Delta T = 0.01$ sec Activate Synchronisation
 Device B 4 samples in $\Delta T = 0.01$ sec

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Activated SYNC Function

Fig. 26 Synchronised data acquisitions

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
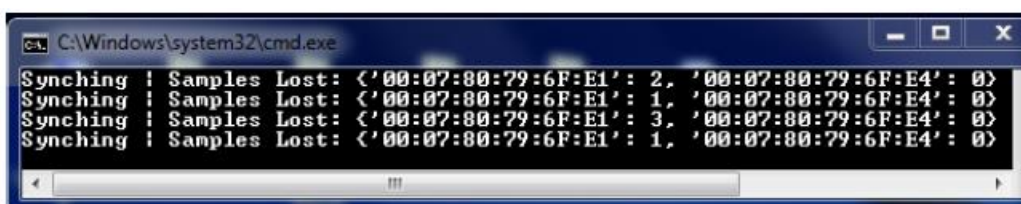
Fig. 27 Open Signals log shown in Windows' command program (CMD.exe)

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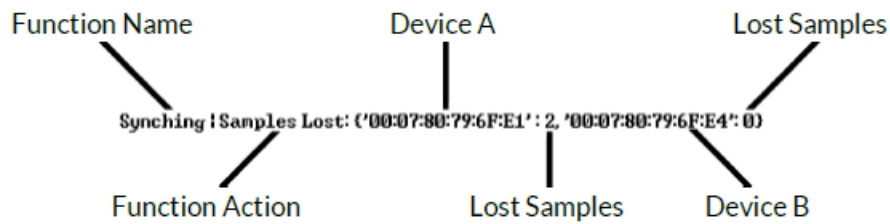- Download the Python API package available at http://downloads.plux.info/PLUX_API_Python2.7_v1.1.zip
- Extract the files from the package to a working folder
- Open the readme.rtf file on the working folder and follow the instructions

9 – Troubleshooting

9.1 – Error List

[Coming Soon]

Error Number: 00000 Error description: [DESCRIPTION]

[SOLUTION]

9.2 – Bluetooth

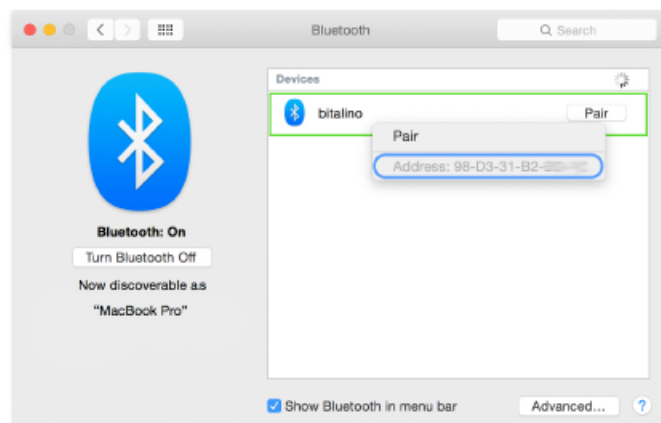
9.2.1 – Which one is my device?

Using several Flux devices at a time, for example using many Bitalinos in a classroom, may be confusing when trying to detect the right device for the right computer. Comparing the device's MAC address to the one shown on your computer will help finding the correct device.

How to show the MAC Address:


Mac OSX – Yosemite and newer

Open the Bluetooth device manager and search for available devices. Right-click on one of the devices to see its MAC address

Mac OSX – Mavericks and older

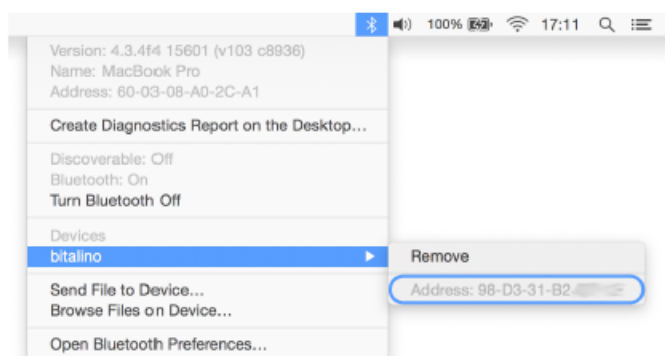Hold the alt key on your keyboard and click on the Bluetooth symbol in the menu bar. A list of the connected devices will appear. Click on your device and check the MAC address on your device with the MAC address on your display.

Windows XP/Vista/7/8

Click the "Start" button and go to the "Control Panel". Select the "Hardware" tab and open the device's properties (right click > properties) and check the shown MAC address with your device's MAC address.

Note: Unfortunately, it is not able to see the device's MAC address before establishing a connection with it in Windows operating systems.

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