

Electrocardiography (ECG) Sensor Data Sheet

ECG 011020

SPECIFICATIONS

- > **Gain:** 1100
- > **Range:** $\pm 1.5\text{mV}$ (with $V_{CC} = 3.3\text{V}$)
- > **Bandwidth:** 0.5-40Hz
- > **Consumption:** $\sim 0.17\text{mA}$
- > **Input Voltage Range:** 2.0-3.5V
- > **Input Impedance:** 7.5GOhm
- > **CMRR:** 86dB

FEATURES

- > Bipolar differential measurement
- > Pre-conditioned analog output
- > High signal-to-noise ratio
- > Small form factor
- > Raw data output
- > Easy-to-use
- > "On-the-person" and "off-the-person" use

APPLICATIONS

- > Heart rate & heart rate variability
- > Human-Computer Interaction
- > Biometrics
- > Affective computing
- > Physiology studies
- > Psychophysiology
- > Biofeedback
- > Biomedical devices prototyping

GENERAL DESCRIPTION

Heartbeats are triggered by bioelectrical signals of very low amplitude generated by a special set of cells in the heart (the SA node). Electrocardiography (ECG) enables the translation of these electrical signals into numerical values, enabling them to be used in a wide array of applications. Our sensors allow data acquisition not only at the chest ("on-the-person"), but also at the hand palms ("off-the-person"), and works both with pre-gelled and most types of dry electrodes. The bipolar configuration is ideal for low noise data acquisition.

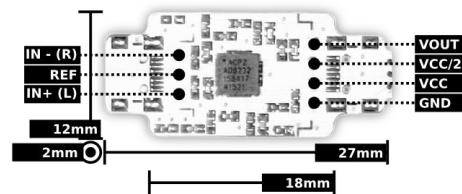


Fig. 1. Pin-out and physical dimensions.

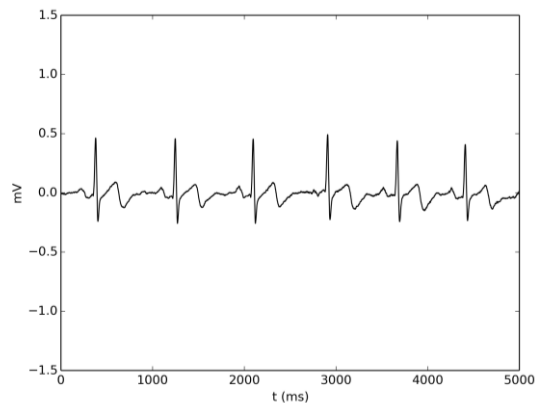


Fig. 2. Typical raw ECG data (acquired with BITalino (r)evolution) using an Einthoven triangle configuration.



Fig. 3. Example of a 1-lead placement with IN+ & IN- on the collarbones and REF on the iliac crest.

bitalino

PLUX – Wireless Biosignals, S.A.
Av. 5 de Outubro, n. 70 – 2.
1050-059 Lisbon, Portugal
bitalino@plux.info
<http://bitalino.com/>

REV B

© 2020 PLUX

This information is provided "as is," and we make no express or implied warranties whatsoever with respect to functionality, operability, use, fitness for a particular purpose, or infringement of rights. We expressly disclaim any liability whatsoever for any direct, indirect, consequential, incidental or special damages, including, without limitation, lost revenues, lost profits, losses resulting from business interruption or loss of data, regardless of the form of action or legal theory under which the liability may be asserted, even if advised of the possibility of such damages.



BEWARE: DIRECT OR INDIRECT COUPLING TO THE MAINS MAY RESULT IN SHOCKING HAZARD



Electrocardiography (ECG) Sensor Data Sheet

TRANSFER FUNCTION

[-1.5mV, 1.5mV]

$$ECG(V) = \frac{\left(\frac{ADC}{2^n} - \frac{1}{2}\right) \times VCC}{G_{ECG}}$$

$$ECG(mV) = ECG(V) \times 1000$$

$VCC = 3.3V$ (operating voltage)

$G_{ECG} = 1100$ (sensor gain)

$ECG(V)$ – ECG value in Volt (V)

$ECG(mV)$ – ECG value in millivolt (mV)

ADC – Value sampled from the channel

n – Number of bits of the channel¹

ORDERING GUIDE

Part #	Description
SENS-ECG-NC	Electrocardiography (ECG) sensor without connectors
SENS-ECG-UCE6	Electrocardiography (ECG) sensor with UC-E6 sockets on both sides for seamless plug & play connection to a BITalino (r)evolution Plugged or Core
SENS-ECG-SHER	Electrocardiography (ECG) sensor with a Molex Sherlock 4-pin socket on one side and a Molex Sherlock 3-pin socket on the other for easy power and signal cable connection or pin breakout using PCB wires

¹ The number of bits for each channel depends on the resolution of the Analog-to-Digital Converter (ADC); in BITalino the first four channels are sampled using 10-bit resolution ($n = 10$), while the last two may be sampled using 6-bit ($n = 6$).