

# Electroencephalography (EEG) Sensor Data Sheet

EEG 26012018

## SPECIFICATIONS

- > **Gain:** 40000
- > **Range:**  $\pm 37.5\mu\text{V}$  (with  $V_{CC} = 3\text{V}$ )
- > **Bandwidth:** 0.8-49Hz
- > **Consumption:**  $\sim 3\text{mA}$
- > **Input Impedance:**  $> 100\text{G}\Omega$
- > **CMRR:** 100dB

## FEATURES

- > Single-channel differential sensor
- > Discrete elastic head band
- > Pre-conditioned analog output
- > High signal-to-noise ratio
- > Shielded miniaturized cables
- > Medical-grade raw data output
- > Ready-to-use form factor

## APPLICATIONS

- > Evoked potentials analysis
- > Neurofeedback
- > Sleep studies
- > Human-Computer Interaction
- > Neurophysiology studies
- > Psychophysiology

## GENERAL DESCRIPTION

Our electroencephalography (EEG) sensor has been especially designed for both classic and localized EEG measurement. When a cap is too intrusive, only a limited number of channels are needed, or you'd like to synchronously record EEG and non-EEG biosignals, this is the perfect solution. The bipolar configuration, with two measurement electrodes detects the electrical potentials in the specific scalp region with respect to a reference electrode, which should be placed in a region of low muscular activity. The resulting signal is the amplified difference between these two signals, eliminating the common unwanted signals detected by the surfaces. Its convenient form factor enables a discrete placement in regions such as the forehead, occipital, and others. Examples:

<http://bit.ly/1E7VenV>  
<http://bit.ly/1PEskAZ>

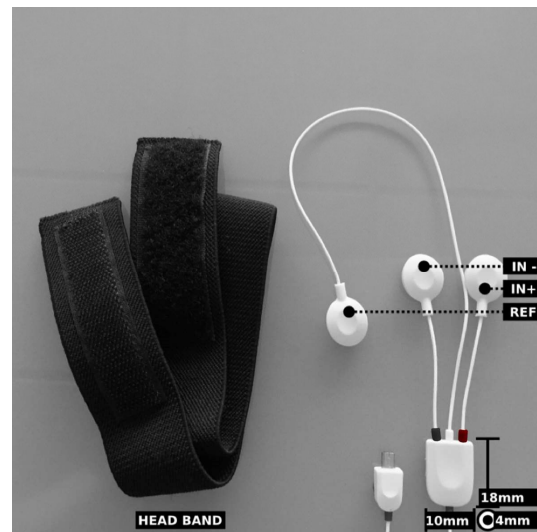


Fig. 1. The sensor is provided with a convenient elastic head band to help secure the electrodes in place.

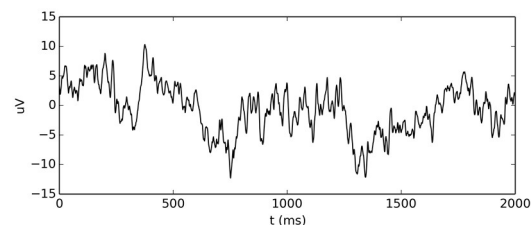


Fig. 2. Typical raw EEG data (acquired with biosignals).

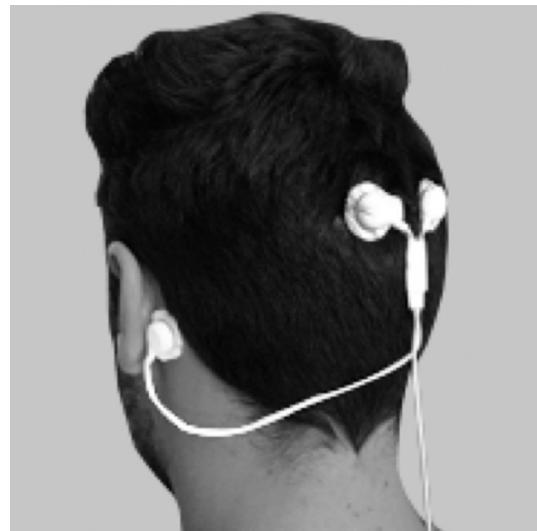


Fig. 3. Example sensor placement for localized EEG.

**biosignalsplux**  
wearable body sensing platForm

PLUX – Wireless Biosignals, S.A.  
Av. 5 de Outubro, n. 70 – 8.  
1050-059 Lisbon, Portugal  
plux@plux.info  
<http://biosignalsplux.com/>

REV A

© 2015 PLUX

This information is provided "as is," and we make no express or implied warranties whatsoever with respect to functionality, operability, use, fitness for a particular purpose, or infringement of rights. We expressly disclaim any liability whatsoever for any direct, indirect, consequential, incidental or special damages, including, without limitation, lost revenues, lost profits, losses resulting from business interruption or loss of data, regardless of the form of action or legal theory under which the liability may be asserted, even if advised of the possibility of such damages.

# Electroencephalography (EEG) Sensor Data Sheet


## PLACEMENT RECOMMENDATIONS



reference electrode placement



sensor electrodes must be 1cm-3cm apart

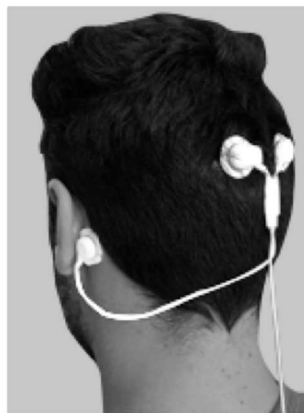

place the sensor on the electrodes




place the elastic band covering the snaps

## USAGE RECOMMENDATIONS

### with noise influence


EEG signal acquisition must be performed in a low electromagnetic noise environment.

A room without power supplies and with the lights off is an appropriate environment to perform the signal acquisitions.

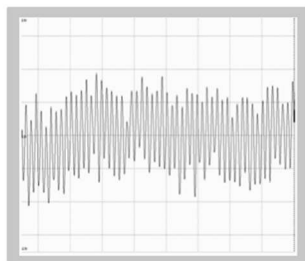


signal acquired in a noisy environment

### our results in a noise controlled room:

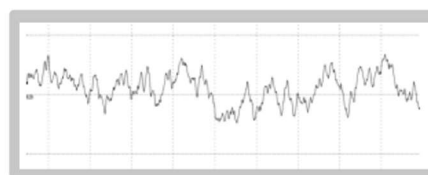

eyes open

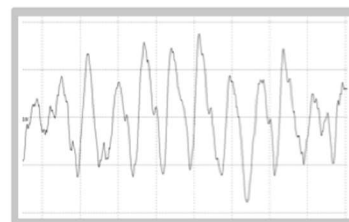

eyes closed

### noise input

The electromagnetic noise enters through the snaps and the cable which connects the sensor to the snaps.

This is a result of not using shielding structures.

# Electroencephalography (EEG) Sensor Data Sheet

---

## TRANSFER FUNCTION

$[-37.5\mu V, 37.5\mu V]$

$$EEG(V) = \frac{\left(\frac{ADC}{2^n} - \frac{1}{2}\right) \cdot VCC}{G_{EEG}}$$

$$EEG(\mu V) = EEG(V) \cdot 1 \times 10^6$$

$VCC = 3V$  (operating voltage)

$G_{EEG} = 40000$  (sensor gain)

$EEG(V)$  – EEG value in Volt (V)

$EEG(\mu V)$  – EEG value in microvolt ( $\mu V$ )

$ADC$  – Value sampled from the channel

$n$  – Number of bits of the channel<sup>1</sup>

## PHYSICAL CHARACTERISTICS

> **W1 x L1 x H1:** 1.0x1.8x0.4cm

> **W2 x L2 x H2:** 1.5x2.3x0.4cm


> **A1:** 105.0±0.5cm

> **A2:** 2.5±0.5cm

> **A3:** 10.0±0.5cm

> **D:** 0.4cm

> **S:** White, Black, Blue, Green, Red, Yellow, Gray, or Brown

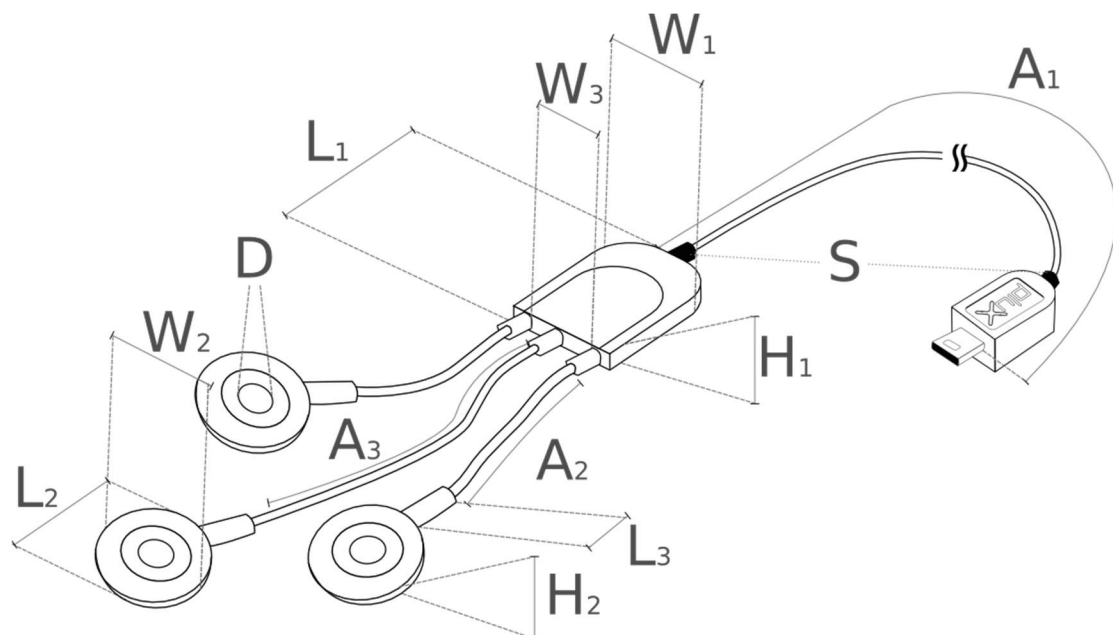

<sup>1</sup> The number of bits for each channel depends on the resolution of the Analog-to-Digital Converter (ADC); in biosignalsplux the default is 16-bit resolution ( $n = 16$ ), although 12-bit ( $n = 12$ ) and 8-bit ( $n = 8$ ) may also be found.

# Electroencephalography (EEG) Sensor Data Sheet

---

## **ORDERING GUIDE**

| <b>Reference</b> | <b>Package Description</b>   |
|------------------|--|
| EEG1             | Electroencephalography (EEG) sensor with standard physical characteristics and a random cable sleeve color   |
| EEG1-A1-A2-A3-S  | Electroencephalography (EEG) sensor built with custom lengths A1, A2 and/or A3 (all in cm), and custom sleeve color S; for standard physical characteristics in A1, A2, A3, or S use 0.<br><br>Examples:<br>> EEG1-200-0-0-0: Otherwise all-standard EEG sensor except for a 200cm cable A1<br>> EEG1-0-0-0-Yellow: Otherwise all-standard EEG sensor except for a yellow cable sleeve<br>> EEG1-50-10-10-Red: Fully custom EEG sensor with a 50cm cable A1, 10cm electrode cables A2 & A3, and a red cable sleeve |