

Light (LUX) Sensor Data Sheet

LUX 090616

SPECIFICATIONS

- > **Range:** 360-970nm (with VCC = 3.3V)
- > **Consumption:** ~50 μ A
- > **Input Voltage Range:** 2.0-6.0V
- > **Angle of Half Sensitivity:** $\pm 60^\circ$

FEATURES

- > Adapted to human eye responsiveness
- > Pre-conditioned analog output
- > High signal-to-noise ratio
- > Small form factor
- > Raw data output
- > Easy-to-use

APPLICATIONS

- > Synchronization with a computer screen
- > Optical marker detector
- > Ambient light monitoring

GENERAL DESCRIPTION

Light (LUX) sensors are typically used for ambient light measurement. However, a common need when working with biosignals is the synchronization of the recorded data with specific light sources (e.g. a computer screen for visual evoked potentials). If applied to the computer screen, our LUX sensor can be used to detect chromatic changes in the stimuli, hence providing a synchronization source. The LUX sensor can also be useful for optical synchronization with third-party devices (provided that such device can trigger an LED), in applications where it is important to have electrical decoupling between devices.

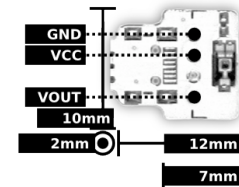


Fig. 1. Pin-out and physical dimensions.

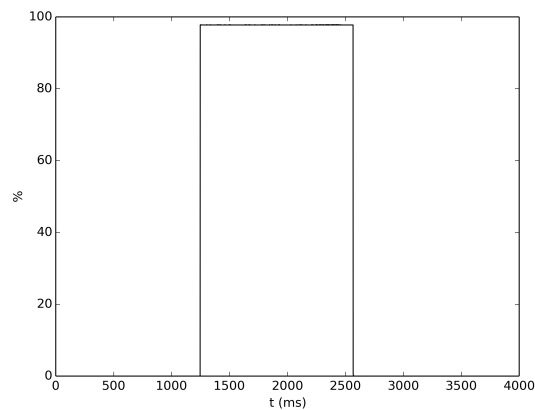


Fig. 2. Typical raw LUX response to a synchronization light source (acquired with BITalino (r)evolution).

bitalino

PLUX – Wireless Biosignals, S.A.
Av. 5 de Outubro, n. 70 – 8.
1050-059 Lisbon, Portugal
bitalino@plux.info
<http://bitalino.com/>

REV A

© 2016 PLUX 

This information is provided "as is," and we make no express or implied warranties whatsoever with respect to functionality, operability, use, fitness for a particular purpose, or infringement of rights. We expressly disclaim any liability whatsoever for any direct, indirect, consequential, incidental or special damages, including, without limitation, lost revenues, lost profits, losses resulting from business interruption or loss of data, regardless of the form of action or legal theory under which the liability may be asserted, even if advised of the possibility of such damages.



BEWARE: DIRECT OR INDIRECT COUPLING TO THE MAINS MAY RESULT IN SHOCKING HAZARD



Light (LUX) Sensor Data Sheet

TRANSFER FUNCTION

[0%, 100%]

$$LUX(\%) = \frac{ADC}{2^n} \cdot 100\%$$

LUX(%) – LUX value in percentage (%)

ADC – Value sampled from the channel

n – Number of bits of the channel¹

ORDERING GUIDE

Part #	Description
SENS-LUX-NC	Light (LUX) sensor without connectors
SENS-LUX-UCE6	Light (LUX) sensor with UC-E6 socket for seamless plug & play connection to a BITalino (r)evolution Plugged or Core
SENS-LUX-SHER4	Light (LUX) sensor with a Molex Sherlock 4-pin socket for easy power and signal cable connection or pin breakout using PCB wires

¹ The number of bits for each channel depends on the resolution of the Analog-to-Digital Converter (ADC); in BITalino the first four channels are sampled using 10-bit resolution ($n = 10$), while the last two may be sampled using 6-bit ($n = 6$).